



PARAFFIN TRIMMER

PARAFFIN TRIMMER BHM1Y1

PARAFFIN TRIMMER



The grooved heating surface channels melted wax to a removable collection tray. The collection tray connects magnetically to the base of the main unit. Disposable plastic liners are included to insert into the collection tray for easy disposal of runoff wax. This unit now includes a digital control panel for personalized settings.

1. Remove the wax trimmer from its packaging and place it on a stable workstation. We recommend that you retain the carton and packaging.
2. The wax trimmer comes with a detachable AC power cord. Connect the cord to the port located on the back of the unit. When ready to use, plug the cord into a power source.
3. Attach the magnetic paraffin collection tray at the bottom of the unit and insert a plastic

SPECIFICATIONS

Model	BHM1Y1
Old Model	BHTP-605
Power Switch & Display	Flip the Power Switch to ON, digital display shows current temperature
Temperature Setting	SET Key to switch to temperature mode, adjust with side buttons, confirm with SET Key
Suggested Operating Temperature	80-90°C (Do not exceed 100°C)
Temperature Control Range	Ambient temperature ~ 110°C
Heating Indicator	Red LED Light indicates heating
Heating Time	Up to 10 minutes to reach desired temperature
Safety Instructions	Do not allow non-cassette items to touch surface, keep sleeves/clothing/jewelry away
Cassette Handling	Cassettes must be flat, pull downward to remove excess paraffin
Net Weight	3 kg
Instrument Dimensions	225*210*140 mm
Package Info	
--Packing	Wooden Case
--Gross Weight	4 kg
--Package Dimensions	370*210*260 mm
Alt Name	Paraffin Trimmer

FEATURES

The grooved heating surface channels melted wax to a removable collection tray. The collection tray connects magnetically to the base of the main unit. Disposable plastic liners are included to insert into the collection tray for easy disposal of runoff wax. This unit now includes a digital control panel for personalized settings.

1. Remove the wax trimmer from its packaging and place it on a stable workstation. We recommend that you retain the carton and packaging.
2. The wax trimmer comes with a detachable AC power cord. Connect the cord to the port located on the back of the unit. When ready to use, plug the cord into a power source.
3. Attach the magnetic paraffin collection tray at the bottom of the unit and insert a plastic collection tray insert. The brim of the plastic liner can be folded down along the side of the tray that sits against the unit. The magnet will hold the collection tray in place.
4. Flip the power switch to the ON position to begin use and adjust temperature. (See Operation section for instructions).



Biolab Scientific Ltd.

Trillium Executive Center, East Tower, 675 Cochrane Dr, Markham, Ontario L3R 0B8, Canada
Email: info@biolabscientific.com | Website: www.biolabscientific.com